

Dissolved Oxygen Polarization Guidelines

More than 30 minutes disconnected from a polarizing device requires 6 hours of polarization.

Less than 30 minutes but greater than 15 minutes disconnected from a polarizing device requires 6 minutes of polarization.

Less than 15 minutes but greater than 5 minutes disconnected from a polarizing device requires 4 minutes of polarization.

Less than 5 minutes disconnected from a polarizing device requires 2 minutes of polarization.

Measurement and Control Products for Science and Industry

19 Thomas, Irvine, California 92618 USA

Phone: (949) 829-5555 Toll-Free: (800) 288-2833 Fax: (949) 829-5560

E-Mail: sales@broadleyjames.com Website: www.broadleyjames.com

