

Don't Worry!

If this flyer doesn't answer your questions, contact our Technical Support:

800-288-2833 (USA & Canada)

Fax: 1-949-829-5560

E-Mail: techsupport@broadleyjames.com

www.biotechcatalog.com




Don't Fret It!

The white crystals you may find on your electrode are formed by the potassium chloride (KCl) in the electrode storage solution. The crystals will not affect the electrode's performance. Rinse the electrode with DI water and proceed as usual.

Don't Sweat It!

The longer the electrode is exposed to high temperatures (above 100°C) the darker it becomes. Likewise, the higher the temperature, the darker the color. This is normal. If the electrode can be calibrated, it's OK to use it.

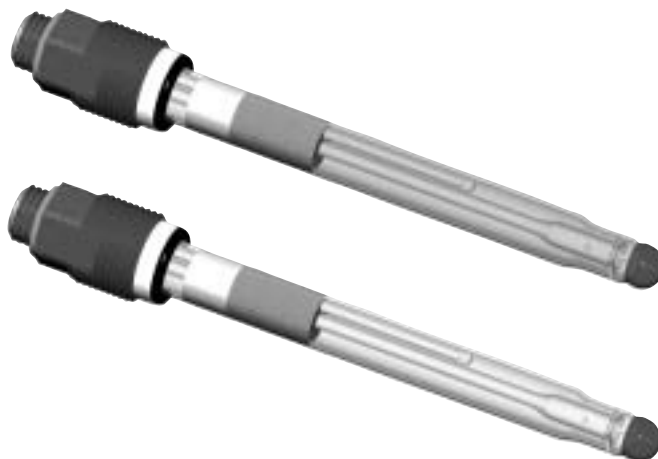

Don't Forget It!

Put the cap on the electrode before autoclaving it. This keeps the connector dry, which insures proper electrode performance.