


# Do you have a datasheet for FermProbe Wetted Materials?

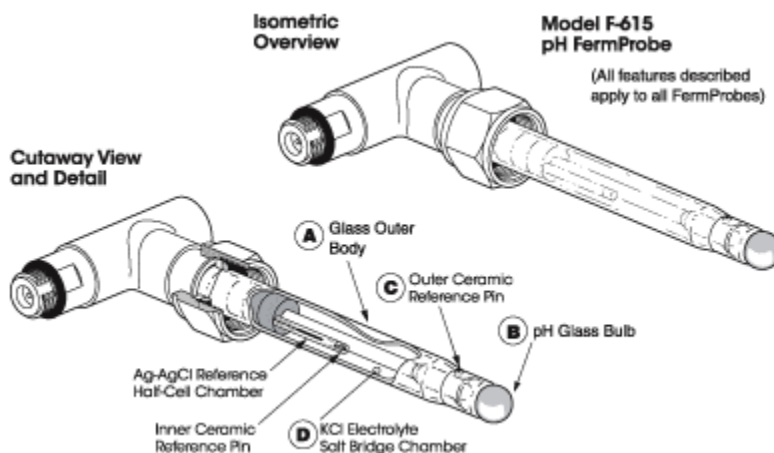

## The FermProbe Double Junction Reference System

All FermProbes have two built-in electrolyte chambers that act to protect and isolate the sensitive inner AgCl reference half-cell. This "double junction", dual-chamber design effectively prevents the most common failure modes of pH electrodes in biopharmaceutical applications.

Only the inner, smaller chamber contains silver ions in the electrolyte. The larger chamber is free of silver ions. This design prevents silver ions from coming into contact with proteins or sulfide ions in the sample media. Reactions between proteins or sulfide ions and the silver ions will cause the formation of substances that will clog the outer ceramic reference pin junction.

All pH and ORP FermProbes have this double junction reference system as a standard feature to insure maximum service life over the widest range of operating conditions.

## Wetted Material List for all FermProbes

- (A) Glass Outer Body: Potash soda lead glass
- (B) Glass Bulb: Silica glass made with small percentages of alkali elements. The exact formulation is proprietary and confidential.
- (C) Porous Ceramic Liquid Junction: Silica plus complex mineral compounds fired at high temperature to form an inert, porous ceramic matrix. The exact formulation is considered proprietary and confidential.
- (D) Internal Electrolyte Gel:

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit [www.broadley-james.com/trademarks](http://www.broadley-james.com/trademarks) for trademark information.

TMP-BF-102101

# Do you have a datasheet for FermProbe Wetted Materials?

1. Potassium Chloride, 3.8 M
2. Cellulose thickening agent, USP Grade

(E) All wetted materials are compliant with applicable FDA regulations.

---

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit [www.broadley-james.com/trademarks](http://www.broadley-james.com/trademarks) for trademark information.

TMP-BF-102101



#### North America and Pacific

Email: [sales@broadleyjames.com](mailto:sales@broadleyjames.com)  
Web: [www.broadleyjames.com](http://www.broadleyjames.com)

19 Thomas, Irvine CA, 92618 USA

Phone: 949-829-5555  
Tollfree: 800-288-2833  
Fax: 949-829-5560

#### United Kingdom and EU Countries

Email: [sales@broadleyjames.co.uk](mailto:sales@broadleyjames.co.uk) Phone: +44 (0)1525 862518  
Web: [www.broadleyjames.eu](http://www.broadleyjames.eu) Fax: +44 (0)1525 862811

Wrest Park, Silsoe Beds MK45 4HS, UK