


How do I change a membrane cartridge on a 19mm and 25mm Probe?

Step 1

While holding the membrane cartridge upright (with opening up) in one hand, carefully pour the electrolyte from the bottle into the cartridge until approximately 75% full. Remove any entrapped air bubbles by tapping on the side of the membrane housing. Try not to spill any of the electrolyte and be certain to wipe off any excess electrolyte with a clean tissue from the outside surfaces of the cartridge.

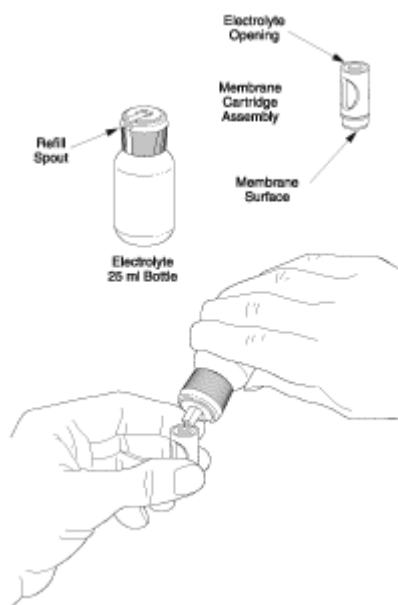

Caution: The use of protective gloves and eyewear is recommended throughout the membrane cartridge refill and installation procedures.

Step 2


In one hand, grip the membrane cartridge, which has been properly filled with electrolyte, by the thumb and forefinger slots while holding the sensor body with the other hand as shown.

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit www.broadley-james.com/trademarks for trademark information.


TMP-BF-102101

How do I change a membrane cartridge on a 19mm and 25mm Probe?

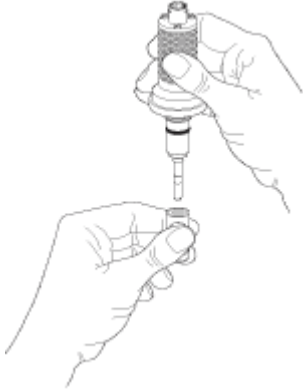

Step 3

While compressing the membrane cartridge with the thumb and forefinger, slide the cartridge over the anode / cathode assembly until the cartridge seats against the conical gasket. Release the pressure from the thumb and forefinger. Rinse lower portion of this subassembly with DI water to remove any excess electrolyte and blot dry with a clean tissue.

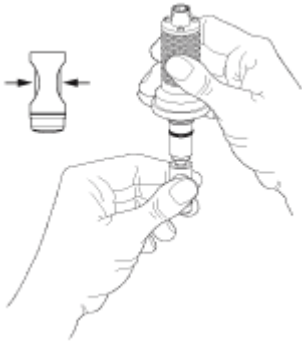

Step 4

The sensor body subassembly with the membrane cartridge in place is now ready for the installation of the stainless steel sensor sleeve.

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit www.broadley-james.com/trademarks for trademark information.


TMP-BF-102101

How do I change a membrane cartridge on a 19mm and 25mm Probe?



Step 5

Slide the stainless steel sensor sleeve over the membrane cartridge, being careful not to nick the o-rings. Be certain that the mating threads of the sleeve and sensor body are free of any dirt or debris.

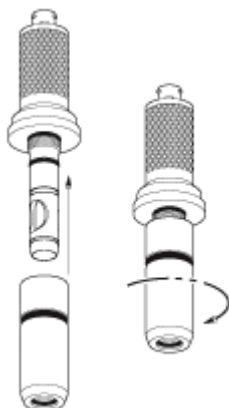

Step 6


Thread the stainless steel sensor sleeve onto the sensor body in the direction shown by the arrow. Be certain that the sleeve is flush with the sensor body so that no o-rings are visible. Do not overtighten. Hand tight is all that is necessary.

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit www.broadley-james.com/trademarks for trademark information.

TMP-BF-102101

How do I change a membrane cartridge on a 19mm and 25mm Probe?

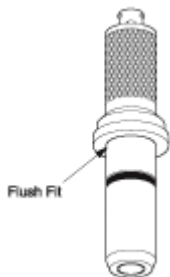

Note: Carefully inspect all o-rings and the conical gasket for any physical damage or excessive wear. Replace these items as needed or if there is any doubt as to their condition.

The contents of this publication are presented for information purposes only, and while every effort has been made to ensure their accuracy, they are not to be construed as warranties or guarantees, express or implied, regarding products or services described herein or in their use or applicability. All sales are governed by our terms and conditions, which are available on request. We reserve the right to modify or improve the design or specification of such products at any time.

© 2022 Broadley-James Corporation. All rights reserved. Visit www.broadley-james.com/trademarks for trademark information.

TMP-BF-102101